

Product Information HA7

Mini Pressure Transmitter HA7

Application/Specified usage

- · Pressure measurement in pipes and vessels
- High Temperature applications up to 150 °C permanent
- Suitable for pressure applications in the Bio-Pharmaceutical and life-science industries

Application examples

- · Ideal for OEM skids such as:
 - Filtration
 - · Chromatography
 - · Blending
 - · CIP/SIP
- Bioreactors
- \cdot Fermenters
- · Pressure vessels
- · Blending and mixing
- · All critical process control

Hygienic design/Process connection

- · Conforms to ASME-BPE 2012 standards
- · CIP-/SIP-cleaning up to 150 °C
- Sensor material and actuating fill are FDA compliant
- · Sensor and product contact surfaces made of stainless steel
- Additional process connections: 3/4", 1", 1¹/₂", 2" Tri-Clamp, Standard & Mini CPM

Features

- · Extremly durable in high temperature applications up to 150 °C permanent
- · Fast response time, less than 200 milliseconds
- · Vacuum-proof
- · Easy to operate
- · Electrical connection with M12 plug-in connector
- · Selectively as relative or absolute measuring sensor available
- Integrated two-wire measurement transmitter 4...20 mA

Options/Accessories

- · Special pressure ranges, customized adjustment ex works
- · Preassembled cable for M12 plug-in connector

Measuring principle of the pressure sensor

This unit utilizes an internal piezoelectric transducer to convert the process measurement into a corresponding mV signal. The mV signal then passes through custom linearization and conditioning circuitry. The resulting signal is an industry standard 4...20 mA. This mA signal is factory set over the specified range of the unit.

With relative (gauge) pressure sensor the back of the transducer is vented to atmosphere, i.e. this sensor measures the gauge pressure and/or vacuum relative to the atmospheric pressure. With absolute pressure sensor the back of the transducer is subject to full vacuum and then permanently sealed, i.e. this sensor measures pressure relative to an absolute vacuum.



PHARMA





ASME BPE 2012

HA7 with Tri-Clamp



PHARMA

Weight

2

Specification		
Pressure ranges	relative [bar] absolute [bar] relative [psi] absolute [psi]	01,0/2,0/3,0/4,0/6,0/7,0/10,0/14,0/20,0/30,0/34,0/35,0/ 50,0/60,0/70,0 -11,0/2,5/3,0/4,0/5,0/6,0/7,0/13,0 02,0/3,0/7,0/14,0/20,0 015/20/30/50/60/100/150/200/300/500/1000 30"Hg0/15/30/60/100/185 020/30/50/60/100/150/200/300
Electrical connection	cable connection supply	M12-plug, 5 pin (1.4305) 1236 V DC
Output	signal burden	420 mA DC, 2 wire loop with non-interrupting circuit verification test points 700 Ω at 24 V
Process	connection	Tri-Clamp 3/4", 1", 1½" and 2" Tri-Clamp Standard and Mini CPM (DIRECTadapt)
Materials	wetted parts housing	stainless steel 1.4404 (316L), R _a ≤ 0.38 µm electropolished stainless steel 1.4305 (304)
Protection class		IP 69 K / NEMA 4X
Accuracy		±0.25 % of full scale
Repeatability		±0.1 % of full scale
Hysteresis		±0.1 % of full scale
Linearity		±0.1 % of full scale
Stability		±0.15 % of calibrated range for 6 months
Temperature ranges	ambient process storage	-1850 °C (0 °F120 °F) -1150 °C (30 °F300 °F) -4065 °C (-40 °F149 °F)
Effect of temperature change		±0.1 psig/10 °F (±7 mbar/5,5 °C) typical
Over-range rating	factor	Minimum of 2 times base range
Response time		< 200 milliseconds
Adjustment	span zero	-50 % of range, except 15 psi / 1 bar rel. and 20 psi / 2 bar abs. cell: -10 % ±10 %

490 g (Tri-Clamp 1.5") 550 g (Tri-Clamp 2") 640 g (CPM) 3



Conventional usage

- · Not suitable for applications in explosive areas.
- · Not suitable for applications in security-relevant equipments (SIL).

PHARMA

Startup

- · Connect the sensor with power supply (12...36 V DC) -> see "Electrical Connection HA7"
- Note: Factory-provided 0...100 % of the complete measuring range are always adjusted to 4...20 mA at the output (for example 0...2 bar at the HA7/057)
- Adjustments of the sensor are made via the potentiometers "ZERO" und "SPAN". Zero point (ZERO) and Span (SPAN) are non-interactive, meaning changing the zero will not change the span.
- · As general maintenance to the unit, a zero check is recommended at approximately 6 month intervals.
- · Other maintenance tasks are not necessary for the pressure sensor HA7.

Electrical Connection / Installation

Although no interaction between zero and span occurs, when making adjustments you should be as close to the top and bottom transmitter range as possible. This will ensure the best possible linearity in the final signal output.

Electrical Connection HA7

Configuration M12-plug

1: + 24 V DC 2: Output 4...20 mA 3: not connected 4: not connected 5: not connected





5

Adjustment of sensor

The standard setting of the HA7 is: 0...100 % of the measurement range (for example 0...2 bar) are equivalent to 4...20 mA at the current output. If it necessary to change these settings for special measurement tasks, perform the following steps:

Adjustment of zero

- Connect ammeter to the current loop (to the terminals TP+ and TP-). If the ammeter displays 4,0 mA no adjustment is necessary.
- In other cases adjust the transmitter "ZERO" until you have the 4,0 mA at the output.
- If you use absolute pressure sensors the expected mA-signal of the zero point has to be calculated in the following way:

mA output = 16 [[Applied pressure – Low end of range] / Transmitter span] + 4 mA

For example:

Applied pressure:	2 bar
Low end of range:	0 bar
Transmitter span:	7 bar
mA output:	8.57 mA

Adjustment of zero



Adjustment of span

Adjustment of span

- · Set the pressure to desired value.
- Connect ammeter to the current loop (to the terminals TP+ and TP-).
- Adjust the transmitter "SPAN" until you have the 20,0 mA you need at the output.
- Factory-provided standard measuring ranges may be turned down to a maximum of 50 % (for example HA7/14REL with measuring range 0...14 bar may be turned down to 0...7 bar)
 Exception:
 15 psi / 1 bar rel. and 20psi / 2 bar abs.: -10 % max.





Cleaning and maintenance

- · Don't use sharp items or aggressive detergents for cleaning.
- · In case of using pressure washers, don't point nozzle directly to electrical connection!

Storage

- No outdoor storage
- · Dry and dust free
- Not exposed to corrosive media
- · Protected against solar radiation
- Avoiding mechanical shock and vibration
- Storage temperature -40...+65 °C
- Relaltive humidity max. 80 %



Transport

- Sensors shall be clean and must not be contaminated with dangerous media!
 - Use suitable transport packaging only to avoid damage of the equipment!

Standards and guidelines



· You have to comply with applicable regulations and directives.

Conformity declaration

Applicable directives:

- EMC Directive 2004/108/EC
- Low Voltage Directive 2006/95/EC
- The CE label confirms compliance of this product with the applicable EC directives.
- · The operator is responsible for ensuring compliance with the directives that are applicable to the overall system.

Disposal

- · This Instrument is not subject to the WEEE directive 2002/96/EG and the respective national laws.
- · Pass the instrument directly on to a specialised recycling company and do not use the municipal collecting points.

Accessories

PVC-cable with M12-connection, 1.4305 (303), IP 69 K, shielded

42117K0025	Molded w/25' cable
42117K0050	Molded w/50' cable
42117K0100	Molded w/100' cable

PVC-cable with M12-connection, 1.4305 (303), IP 69 K, shielded to nut

42117H0025	5-Conductor Molded w/25' cable
42117H0050	5-Conductor Molded w/50' cable
42117H0100	5-Conductor Molded w/100' cable









7

PHARMA

8

Order code

HA7	Compact	pressure t	ransmitte	r				
	Measuren Compatib 246 057 235 192 060 309 061 062 065 066 491 224 068 069 206	nent range le with rar (01 bar) (02 bar) (03 bar) (04 bar) (04 bar) (07 bar) (010 ba (014 ba (014 ba (030 ba (030 ba (035 ba (050 ba (060 ba (070 ba	e bar (0U nge types E)))))) r) (range B r) (range B r) (range B r) (range B r) (range B r) (range B r) (range B	RL) 3 & T ur only) only) only) only) only) only)	iless n	oted	Measu Compa 063 065 066 068 069 071 073 075 077 081 084	rement range psi (0URL) atible with range types G & A unless noted (015 psi) (020 psi) (030 psi) (050 psi) (060 psi) (0100 psi) (0100 psi) (0200 psi) (0200 psi) (0300 psi) (0500 psi) (range G only) (01000 psi) (range G only)
	Measuren Compatib 418 251 286 217 056 356 475 304 495	nent range le with rar (-10 bar (-11 bar (-12.5 b (-13 bar (-14 bar (-15 bar (-16 bar (-17 bar (-17 bar (-113 bar (-17 bar (-113 bar (-113 bar (-113 bar (-113 bar (-113 bar (-113 bar) (-113 bar (-113 bar) (-113 bar (-113 bar) (-113 bar) (-11	e bar (-1l nge types E r) r) r) r) r) r) r) r) r) r) r) r) r)	e (no va ive ute tive onnect (Tri-Cl (Tri-Cl (Tri-Cl (Tri-Cl (Tri-Cl (CPM)) (mini Senso 0	cuum) 5 only ion/fit lamp 1 lamp 1 lamp 2 CPM) or type Stand Diapł 1) tting (/4") ") 1/2") tard hragm m 316L S Finish 3	Measu Compa 025 028 029 031 032 494	rement range psi (30"HgURL) titble with range types C unless noted (30"Hg0) (range M only) (30"Hg30 psig) (30"Hg100 psig) (30"Hg100 psig) (30"Hg185 psig) (30"Hg185 psig)
HA7	7461	B /	004/	0 /	1/	3/	A0 A1 ↓	(QDR & Field wirable connector (FWC) with no cable)

50055 / 2.0 / 2015-02-16 / MU / EU

NEGELE MESSTECHNIK GMBH Raiffeisenweg 7 87743 Egg an der Guenz Phone +49 (0) 83 33 . 92 04 - 0 Fax +49 (0) 83 33 . 92 04 - 49 sales@anderson-negele.com Tech. Support: support@anderson-negele.com Phone +49 (0) 83 33 . 92 04 - 720